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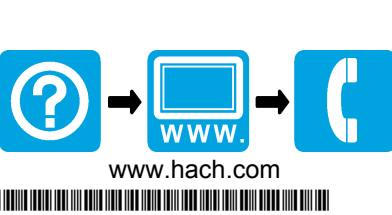
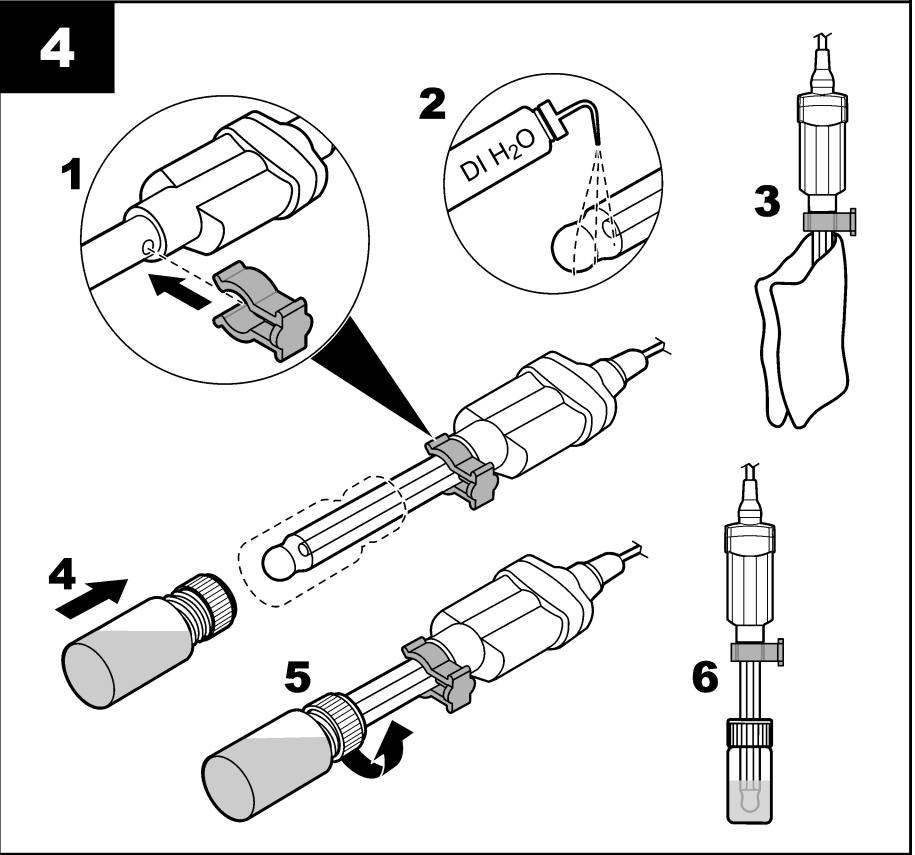
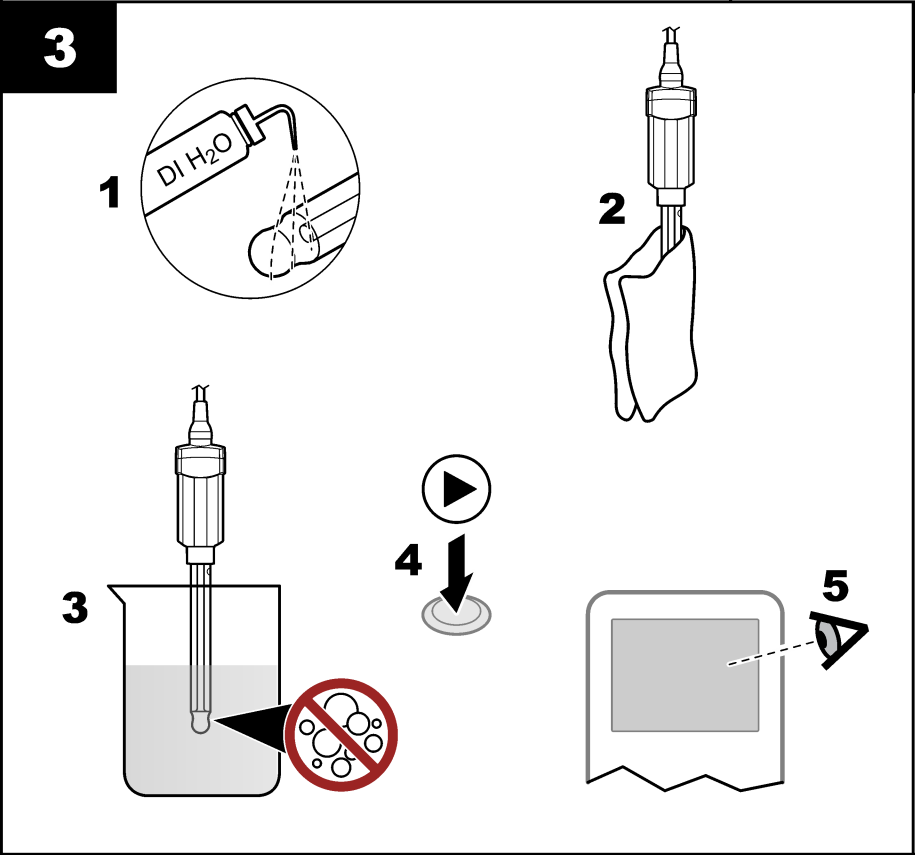
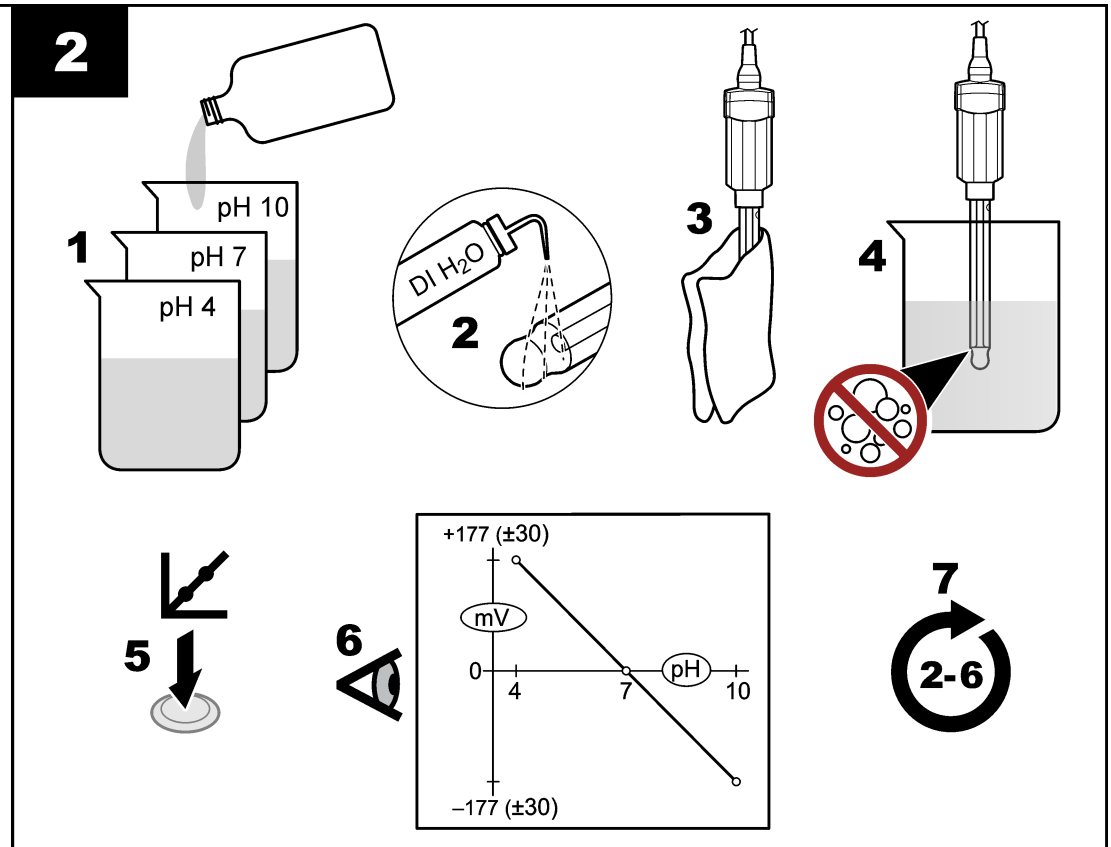
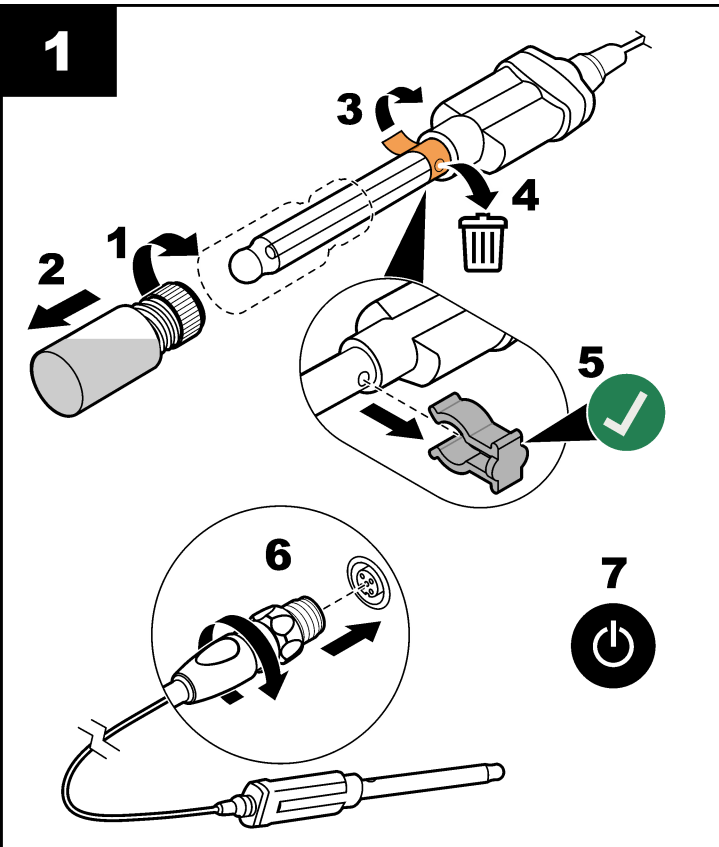
PHC805

PHC805
Introduction

The probe is a pH electrode with refillable electrolyte, a glass body and an integrated temperature sensor. The probe measures pH in general aqueous samples.

Warning
 ⚠ *Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.*

User manual
 A user manual with all instructions is on the manufacturer's website.

Items to collect:
 Compatible meter, pH standard solutions, beakers, DI water, no-lint cloth, filling solution

- 1 Connect the probe**
- 2 Calibration**
Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.
- 3 Measurement**
- 4 Storage**

Technical information

pH range	pH 0 to 14
Slope	-59 mV/pH (85 to 115% at 25 °C)
Temperature	0 to 80 °C (32 to 176 °F)
Minimum sample depth	18 mm (0.71 in.)
Filling solution	3 M KCl
Storage solution	3 M KCl

Maintenance and storage
 For maintenance and storage details, refer to the expanded version of this manual on the manufacturer's website.

Warranty
 1 year for manufacturing faults. This warranty does not cover improper use or wear.

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